2000005301

Heating block_102015

IKA®

Ursprungssprache

DE

IKA® Heizblock IKA® Heating block

Technische Information
Technical information

DE
EN

IKA®-Werke, Germany
Reg. No. 004343

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Explication of warning symbols

-  **DANGER** Indicates an (extreme) hazardous situation, which, if not avoided, will result in death, serious injury.
-  **WARNING** Indicates a hazardous situation, which, if not avoided, can result in death, serious injury.
-  **CAUTION** Indicates a potentially hazardous situation, which, if not avoided, can result in injury.
-  **NOTICE** Indicates practices which, if not avoided, can result in equipment damage.
-  **WARNING** Reference to exposure to a hot surface!
-  Hot surface!

Safety instructions

- Read the operating instructions in full before starting up and follow the safety instructions.
- Keep the operating instructions in a place where it can be accessed by everyone.
- Ensure that only trained staff work with the equipment.
- Follow the safety instructions, guidelines, occupational health and safety and accident prevention regulations.
- The heating blocks maximum allowed temperature must not exceed 300 °C.

 **WARNING** During handling the heating blocks. Hot surface can cause risk of burning! Wear gloves when handling the heating blocks!

- Pay attention to the residual heat after processing.
- Never move the heating blocks as long as they are hot.
- Observe the operating instructions of the respective magnetic stirrer.

 **WARNING** Beware of hazards due to:

- flammable materials
- combustible media with a low boiling temperature
- glass breakage
- incorrect container size
- unsafe condition of container.

 **NOTICE** The round bottom flask must be firmly clamped when it is used with the heating block.

Proper use


The **IKA® Heating block** has been designed for **IKA®** magnetic stirrers with heating function with a maximum working plate diameter 135 mm. The **Heating block** allows syntheses to be carried out in small reaction vessels or round flasks. This system ensures optimal heat transfer from the heating plate into the medium.


Note: Please don't use silicone oil or water as heat transfer medium between carrier and inlay since the inlay or the vessel might rotate inside the carrier block. The anodizing color of the heating blocks could change permanently when temperature is higher than 180 °C or with aging. This does not affect the heating function of the block.

Glassware checklist

- Round bottom flasks or small reaction vessels used with heating blocks/inserts must be free of cracks. Etched or severely scratched glassware should not be used for heating with blocks or inlays.
- We recommend to use round bottom flask that comply with DIN ISO 1773:1999-05 dimension specification. Flask diameters which do not comply with ISO 1773:1999-05 should not exceed the following values:

Flask capacity	Max. diameter
10 ml	33.5 mm
25 ml	43.5 mm
50 ml	52.5 mm
100 ml	66 mm
250 ml	87.5 mm
500 ml	107.5 mm
1000 ml	134.5 mm
2000 ml	169.5 mm
3000 ml	198.5 mm
5000 ml	230.5 mm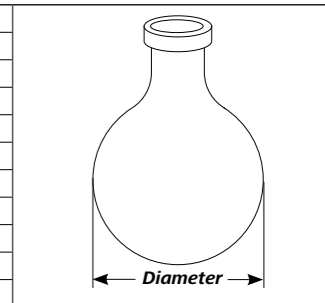

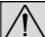 **NOTICE** Use of glassware which exceed the maximum allowed diameter can cause breakage of the glassware!

Cleaning

 **WARNING** Pay attention to the residual heat prior to cleaning. Never clean the devices when they are still hot!

Use only cleaning agents which have been approved by **IKA®** to clean the device.

Dirt	Cleaning agent
Dye	Isopropyl alcohol
Construction material	Water containing tenside / isopropyl alcohol
Cosmetics	Water containing tenside / isopropyl alcohol
Foodstuffs	Water containing tenside
Fuel	Water containing tenside

For materials which are not listed, please request information from **IKA®** application support. Wear protective gloves while cleaning the device. If a different cleaning or decontamination method than the method defined by **IKA®** is planned, the user must ascertain with **IKA®** that this method does not damage the device.


Note: Please don't clean the devices in a lab dishwasher! Or else, the coating may peel off, but has not affect to the functionality of the device.

Warranty

In accordance with **IKA®** warranty conditions, the warranty period is 24 months. For claims under the warranty please contact your local dealer. You may also send the machine direct to our factory, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs. The warranty does not cover worn out parts, nor does it apply to faults resulting from improper use, insufficient care or maintenance not carried out in accordance with the instructions in this operating manual.

IKA®-Werke GmbH & Co. KG
 Janke & Kunkel-Str. 10
 D-79219 Staufen
 Tel. +49 7633 831-0
 Fax +49 7633 831-98
 sales@ika.de

www.ika.com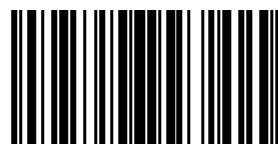

4441200d