2000009407

H 1000...8000_012016

IKA®

IKA® H 1000
IKA® H 1500
IKA® H 3000
IKA® H 5000
IKA® H 8000

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Explication of warning symbols

-  **DANGER** Indicates an (extreme) hazardous situation, which, if not avoided, will result in death, serious injury.
-  **WARNING** Indicates a hazardous situation, which, if not avoided, can result in death, serious injury.
-  **CAUTION** Indicates a potentially hazardous situation, which, if not avoided, can result in injury.
-  **NOTICE** Indicates practices which, if not avoided, can result in equipment damage.
-  **WARNING** Reference to exposure to a hot surface!

Safety instructions

- Read the operating instructions in full before starting up and follow the safety instructions.
- Keep the operating instructions in a place where it can be accessed by everyone.
- Ensure that only trained staff work with the equipment.
- Follow the safety instructions, guidelines, occupational health and safety and accident prevention regulations.

-  **WARNING** During heating, the handles can be very hot.
-  **CAUTION** Never use the beakers with damaged handles.
-  **WARNING** During experiment, exercise caution when handling the beakers. Only touch the beakers at the handles.
 - Pay attention to the residual heat after processing.
 - Never move the beakers as long as they are hot.
 - Observe the operating instructions of the respective magnetic stirrer.
 - The beaker must be properly placed on the heating plate and must not come off by itself. The gravity centre of beaker must lie in the heating plate.
-  **WARNING** Beware of hazards due to:
 - flammable materials
 - combustible media with a low boiling temperature
 - glass breakage
 - incorrect container size
 - unsafe condition of container.

Proper use

The beakers can be used for matching with **IKA**® magnetic stirrers with heating or heating plates. **H 3000**, **H 5000** and **H 8000** beaker has round deepening area for fitting heating plates having a diameter of 135 mm. The stainless steel beakers ensure good heat transfer and quick heating-up of tempering medium. The handles protect you from burning caused by hot medium.

Cleaning

 **WARNING** Pay attention to the residual heat prior to cleaning. Never clean the devices when they are still hot!

Use only cleaning agents which have been approved by **IKA**® to clean the device.

Dirt	Cleaning agent
Dye	Isopropyl alcohol
Construction material	Water containing tenside / isopropyl alcohol
Cosmetics	Water containing tenside / isopropyl alcohol
Foodstuffs	Water containing tenside
Fuel	Water containing tenside

For materials which are not listed, please request information from **IKA**® application support. Wear protective gloves while cleaning the device. If a different cleaning or decontamination method than the method defined by **IKA**® is planned, the user must ascertain with **IKA**® that this method does not damage the device.

Technical data

		H 1000	H 1500	H 3000	H 5000	H 8000
Inner diameter	mm	160	140	180	220	265
Height	mm	79	160	119	142	162
Volume	ml	1000	1500	3000	5000	8000
Fit heating plate size	mm	no size limitation			Ø 135	
Material		Stainless steel 1.4301 (AISI 304)				


Subject to technical changes!

Warranty

In accordance with **IKA**® warranty conditions, the warranty period is 24 months. For claims under the warranty please contact your local dealer. You may also send the machine direct to our factory, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs. The warranty does not cover worn out parts, nor does it apply to faults resulting from improper use, insufficient care or maintenance not carried out in accordance with the instructions in this operating manual.

IKA®-Werke GmbH & Co.KG
 Janke & Kunkel-Str. 10
 D-79219 Staufen
 Tel. +49 7633 831-0
 Fax +49 7633 831-98
 sales@ika.de

www.ika.com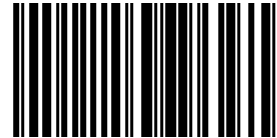

25001879a